


LOOXster®

Enabling technology
for bacterial and fungal nucleic acid detection

- Enrichment of bacterial or fungal DNA
- Reduction of host background DNA
- Improvement of performance and sensitivity of downstream protocols

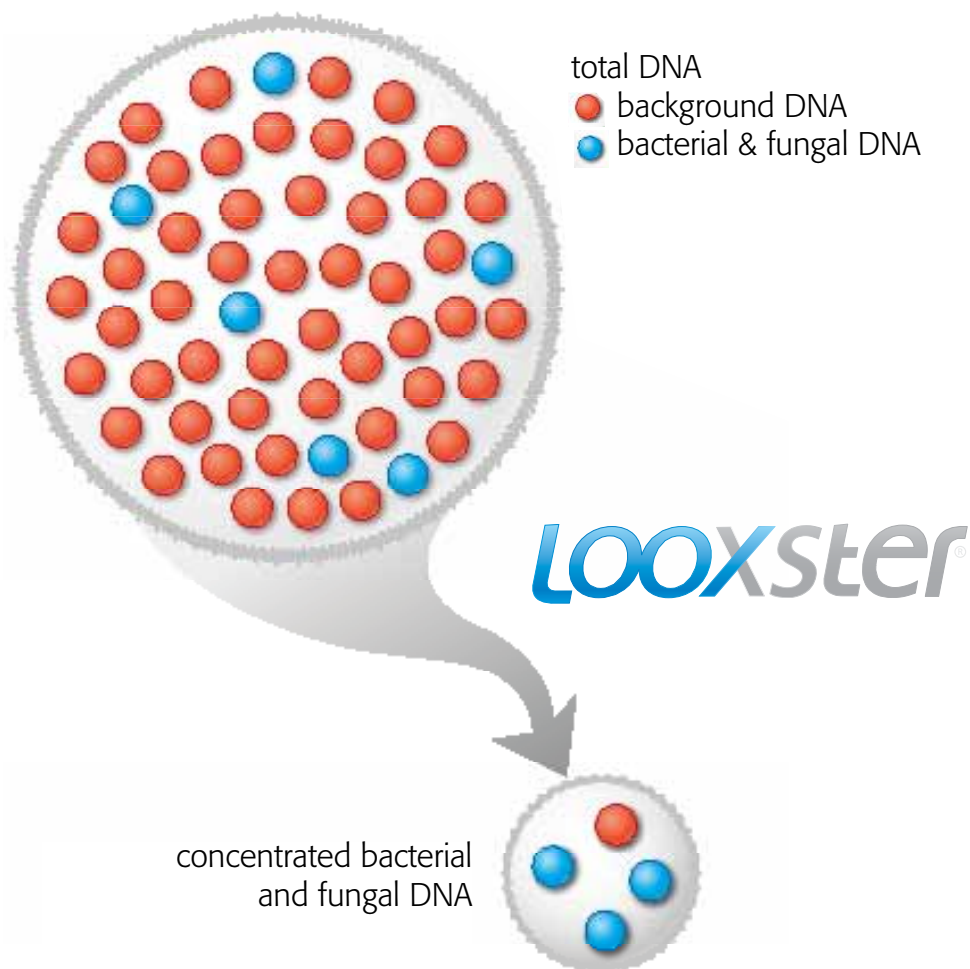

Our value proposition


LOOXSTER® is an enabling technology that makes DNA analyses more sensitive and methods for DNA preparation more effective.

What does LOOXSTER® do?

LOOXSTER® is able to enrich prokaryotic and fungal DNA molecules in samples containing an excess of higher eukaryotic DNA. Due to the reduction of most of the background DNA LOOXSTER® makes large sample sizes accessible to analytical and preparative downstream applications.

How can LOOXSTER® help you?

In qPCR target DNA concentration is crucial for reliable and sensitive analyses. To demonstrate the LOOXSTER®-effect 50 µg of human genomic DNA isolated from an EDTA blood sample was spiked with 10³ copies of *E.coli* DNA and 200 ng of LOOXSTER®-treated and untreated template DNA was used for qPCR. The comparative qPCR shows a ΔCt of 5,23 revealing a 37,5 fold increase in target DNA concentration in the LOOXSTER®-sample compared to the untreated counterpart.

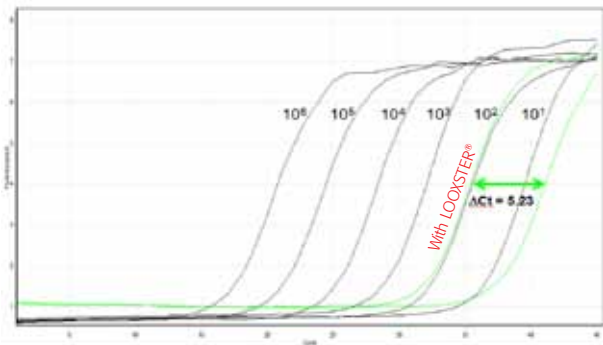

Comparative qPCR

Features


- Enrichment of prokaryotic and fungal DNA in samples containing an excess of higher eukaryotic DNA
- Removal of more than 95 % of higher eukaryotic DNA
- Up to 300 µg of input-DNA
- DNA clean-up included

How does LOOXSTER® work?

LOOXSTER® uses the specific affinity of the CXXC finger protein 1 (CFP1) for non-methylated CpG-dinucleotides^[1]. DNA containing non-methylated CpG-dinucleotides binds efficiently to LOOXSTER® and can therefore be separated from DNA that underwent CpG methylation^[2]. In general, the efficiency of LOOXSTER® DNA enrichment depends on differential DNA G+C contents and DNA methylation rates among species.

What are typical fields of LOOXSTER® application?

PCR and qPCR, array analysis, DNA sequencing, library construction etc. in pathogen detection^[3-7], food and biosafety, biopharma, molecular ecology and microbiome research.

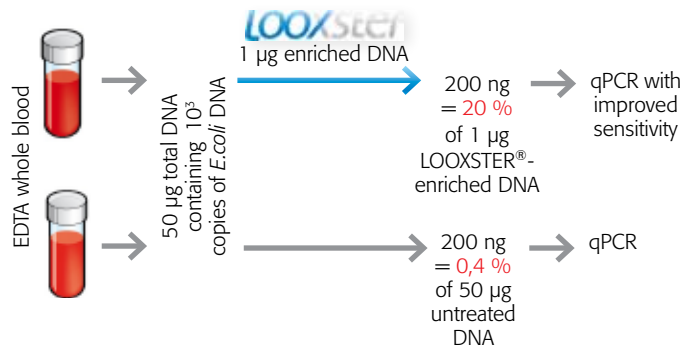

LOOXSTER® principle

Order information

LOOXSTER® Blood & Tissue DNA Kit - KFFLX

Complete sample prep comprising mechanical lysis, DNA extraction and LOOXSTER® enrichment, operated on a KingFisher® FLEX instrument.

| Order number | Quantity |
|----------------|--------------|
| 847-0209500401 | 12 reactions |
| 847-0209500402 | 24 reactions |

LOOXSTER® Enrichment Kit

Manual LOOXSTER® enrichment from up to 300 µg total DNA preparation.

| Order number | Quantity |
|----------------|--------------|
| 847-0209500302 | 10 reactions |

LOOXSTER® is for research use only. LOOXSTER® is a registered trademark.

References

1. Voo K S, Carlone D L, Jacobsen B M, Flodin A, Skalnik D G. Cloning of a mammalian transcriptional activator that binds unmethylated CpG motifs and shares a CXXC domain with DNA methyltransferase, human trithorax, and methyl-CpG binding domain protein 1. *Mol Cell Biol.* 2000;20:2108-2121.
2. Sachse, S., E. Straube, M. Lehmann, M. Bauer, S. Rußwurm, K.-H. Schmidt. 2009. Truncated hCGBP improves NAT-based detection of bacterial species in human samples. *J Clin Microbiol* 47(4):1050-1057.
3. Kriegsmann, J., T. Hopf, D. Jacobs, N. Arens, V. Krenn, R. Schmitt-Wiedhoff, M. Kriegsmann, M. Heisel, C. Biehl, H. Thabe, R.P.H. Schmitz, M. Lehmann, M. Otto. 2009. Molekularpathologische Diagnostik von Gelenkinfektionen. *Orthopade* 38(6):531-538.
4. Sachse, S., M. Lehmann, R.P.H. Schmitz, K. Reinhart, E. Straube. 2008. Multiplex PCR vs. Blood Culture: Towards a New Gold Standard in Sepsis Diagnostics. International Shock Congress, 6th Congress of the International Federation of Shock Societies, 31st Annual Conference on Shock, and 7th International Conference on Complexity in Acute Illness. 28.06.-02.07.2008, Cologne, Medimond Proceedings, Bologna, Italy
5. Wiesinger-Mayr H, Jordana-Lluch E, Martró E, Schoenthaler S, Noehammer C. Establishment of a semi-automated pathogen DNA isolation from whole blood and comparison with commercially available kits. *J Microbiol Methods.* 2011 Jun;85(5):206-13. doi: 10.1016/j.mimet.2011.03.003. Epub 2011 Mar 21.
6. Horz HP, Scheer S, Vianna ME, Conrads G. New methods for selective isolation of bacterial DNA from human clinical specimens. *Anaerobe.* 2010 Feb;16(1):47-53. doi: 10.1016/j.anaerobe.2009.04.009. Epub 2009 May 20.
7. Horz HP, Scheer S, Huenger F, Vianna ME, Conrads G. Selective isolation of bacterial DNA from human clinical specimens. *J Microbiol Methods.* 2008 Jan;72(1):98-102. Epub 2007 Nov 28.