





innuPREP DNA Mini Kit

Protocol 2: DNA extraction from paraffin embedded tissue samples



- Recommended steps before starting
- Heat thermal mixer or water bath (37 °C/ 50 °C/ 90 °C)
 - Prepare Proteinase K, Washing Solution HS and Washing Solution MS according to the instruction

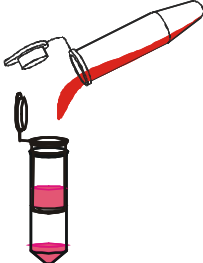

-
1. Starting material
- Place paraffin embedded tissue samples into a 2.0 ml tube



-
2. Remove paraffin
- 
- 
- Add 1 ml Octane or Xylene, vortex
 - Centrifuge: max speed, 5 min
 - Discard supernatant

-
3. Washing (Repeat 2 x)
- 
- 
- Add 1 ml ethanol (96 – 99.8 %)
 - Vortex
 - Centrifuge: max speed, 3 min
 - Remove ethanol (pipetting)

-
4. Evaporate residual ethanol
- Incubation: 10 – 15 min @ 37 °C

-
5. Lysis
- 
- 
- Add 400 µl TLS and 25 µl PK
 - Vortex: 5 sec
 - Incubation: until sample is lysed @ 50 °C
 - Incubation: 60 min @ 90 °C

-
6. Binding of DNA
- 
- 
- Add 400 µl TBS
 - Vortex: 15 sec
 - Add Spin Filter to Receiver Tube
 - Add sample to Spin Filter
 - 11,000 x g (~11,000 rpm): 2 min

-
7. Washing
New Receiver Tube
- 
- 
- Add 500 µl HS
 - 11,000 x g (~11,000 rpm): 1 min
 - Add 750 µl MS
 - 11,000 x g (~11,000 rpm): 1 min

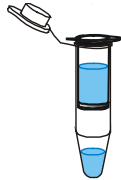
8. Remove Ethanol

New Receiver Tube



- Discard filtrate
- Add Spin Filter to Receiver Tube
- 11,000 x g (~11,000 rpm): 2 min

9. Elution



- Add Spin Filter to Elution Tube
- Add 200 µl Elution Buffer
- Incubation: 1 min @ RT
- 11,000 x g (~11,000 rpm): 1 min

Order No.:	845-KS-1041010	10 reactions
	845-KS-1041050	50 reactions
	845-KS-1041250	250 reactions

This documentation describes the state at the time of publishing. It needs not necessarily agree with future versions. Subject to change!

Expression and further use permitted with indication of source. © 2018 Analytik Jena AG, AJ Innuscreen GmbH

Analytik Jena AG
Konrad-Zuse-Straße 1
07745 Jena · Germany

Phone +49 3641 77 9400
Fax +49 3641 77 767776
info@analytik-jena.com
www.analytik-jena.com

Manufacturer:
AJ Innuscreen GmbH
Robert-Rössle-Straße 10
13125 Berlin