



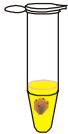

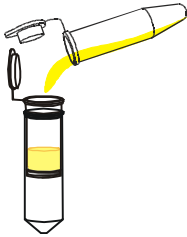




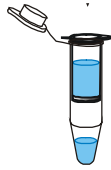
innuPREP DOUBLEpure Kit

Protocol: DNA extraction from agarose gel slices

- Recommended steps before starting
- Heat thermal mixer or water bath (50 °C)
 - Prepare Washing Solution LS according to the instruction

-
1. Starting material
- TAE agarose gels
 - TBE agarose gels
 - Up to 300 mg
 - Up to 300 mg
-
2. Lysis
- 
- Add 650 µl Gel Solubilizer
 - Incubation: 50 °C, 10 min
-
3. Binding of DNA
- Re-use Receiver Tube
- 
- 
- Add 50 µl Binding Optimizer
 - Mix: vortex or pipetting
 - Add Spin Filter to Receiver Tube
 - Add sample to Spin Filter
 - 11,000 x g (~11,000 rpm): 1 min
-
4. Washing
- Re-use Receiver Tube
- 
- 
- Add 700 µl Washing Solution LS
 - 11,000 x g (~11,000 rpm): 1 min
 - Add 700 µl LS
 - 11,000 x g (~11,000 rpm): 1 min
-
5. Remove Ethanol
- Re-use Receiver Tube
- 
- 
- Discard filtrate
 - Add Spin Filter to Receiver Tube
 - Centrifuge: max speed, 2 min
-
- 6.1 Elution (standard elution volume)
- 
- 
- Add Spin Filter to an Elution Tube
 - Add 30–50 µl Elution Buffer
 - Incubation: 1 min @ RT
 - 11,000 x g (~11,000 rpm): 1 min
-

6.2 Elution
(mini elution
volume)



- Add Spin Filter to an Elution Tube
- Add 10–20 µl Elution Buffer
- Incubation: 2 min @ RT
- 11,000 x g (~11,000 rpm): 1 min

Order No.:	845-KS-5050010	10 reactions
	845-KS-5050050	50 reactions
	845-KS-5050250	250 reactions

This documentation describes the state at the time of publishing. It needs not necessarily agree with future versions. Subject to change!

Expression and further use permitted with indication of source. © 2021 Analytik Jena GmbH, AJ Innuscreen GmbH

Analytik Jena GmbH
Konrad-Zuse-Straße 1
07745 Jena
Germany

Phone +49 3641 77 94 00
Fax +49 3641 77 767776
info@analytik-jena.com
www.analytik-jena.com

Manufacturer:
AJ Innuscreen GmbH
Robert-Rössle-Straße 10
13125 Berlin