


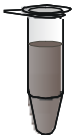






innuPREP Plasmid Mini Kit

Protocol 2: Isolation of plasmid DNA from 5-15 ml bacterial culture

Recommended steps before starting

- Prepare Washing Solution B according to the instruction

1. Starting material	<i>E. coli</i> overnight culture		<ul style="list-style-type: none"> ▪ > 5–15 ml
2. Pellet cells			<ul style="list-style-type: none"> ▪ Centrifuge: max. speed, 8 min ▪ Discard supernatant completely
3. Resuspend cells			<ul style="list-style-type: none"> ▪ Add 550 µl Resuspension Buffer ▪ Vortex ▪ Transfer solution to 2.0 ml tube
4. Lysis Don't vortex!			<ul style="list-style-type: none"> ▪ Add 550 µl Lysis Buffer ▪ Mix: invert tube 6–8 times ▪ Lysis time: ≤ 5 min
5. Neutralization			<ul style="list-style-type: none"> ▪ Add 750 µl Neutralization Buffer ▪ Mix: invert tube 6–8 times ▪ Centrifuge: max speed, 8 min
6. Binding of DNA Re-use Receiver Tube		 	<ul style="list-style-type: none"> ▪ Add Spin Filter to Receiver Tube ▪ Add 850 µl of clarified sample to Spin Filter ▪ 11,000 x g (~11,000 rpm): 1 min ▪ Discard filtrate and load residual sample to Spin Filter ▪ 11,000 x g (~11,000 rpm): 1 min

7. Washing

Re-use Receiver Tube



- Add 650 μ l Washing Solution A
- 11,000 x g (~11,000 rpm): 1 min
- Add 750 μ l Washing Solution B
- 11,000 x g (~11,000 rpm): 1 min

8. Remove Ethanol

Re-use Receiver Tube



- Discard filtrate
- Add Spin Filter to Receiver Tube
- Centrifuge: max speed, 3 min

9. Elution



- Add Spin Filter to a 1.5 ml reaction tube
- Add 100 μ l Elution Buffer P
- Incubation: 3 min @ RT
- 11,000 x g (~11,000 rpm): 1 min

Order No.:	Spin Filter with cap
	845-KS-5041010 10 reactions
	845-KS-5041050 50 reactions
	845-KS-5041250 250 reactions
	845-KS-5041500 500 reactions

This documentation describes the state at the time of publishing. It needs not necessarily agree with future versions. Subject to change!

Expression and further use permitted with indication of source. © 2021 Analytik Jena GmbH, AJ Innuscreen GmbH

Analytik Jena GmbH
Konrad-Zuse-Strasse 1
07745 Jena · Germany

Phone +49 3641 77 70
Fax +49 3641 77 9279
info@analytik-jena.com
www.analytik-jena.com

Manufacturer:
AJ Innuscreen GmbH
Robert-Rössle-Strasse 10
13125 Berlin · Germany